


PS-3215; Wireless Colorimeter Sensor Battery Replacement Instructions

Reference: Battery part number 540-113; Li-Poly 1000 mAH

Recommended Equipment: (PS-3299; Battery Replacement Tool Set)




Torx T8 Driver

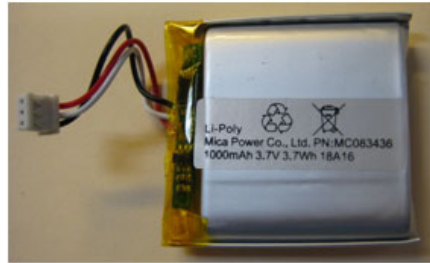

Replacement Battery



CAUTION: Failure to use proper ESD protection could result in damage to the electronics.

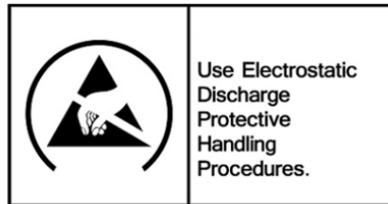

Step 1: Position the Sensor on a surface as shown.

- Peel off the (4) rubber feet using any small flat blade screwdriver.
- Place the rubber feet with the sticky side facing upward on the surface.
- Remove the (4) T8 Screws with a Torx T8 Screwdriver.





Position as Shown



Remove Rubber Feet

Remove Qty=4 Screws

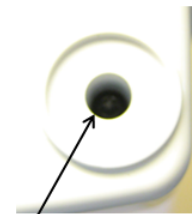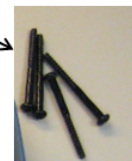


Torx T8 Screws (4)

PASCO Form Number	Description	Date	Page	Revision
013-16110	PS-3215 Battery Replacement Instructions	05-05-22	Page 1 of 3	B

Step 2: While holding both halves together, flip the product over as shown.




Step 3: Lift off the top cover and place it as shown.

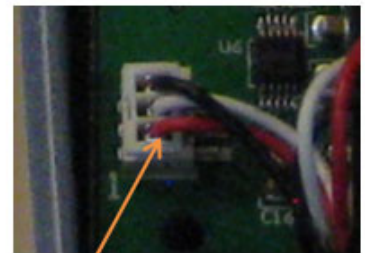

Battery Connector
Note: red wire
location

PASCO Form Number	Description	Date		Revision
013-16110	PS-3215 Battery Replacement Instructions	05-05-22	Page 2 of 3	B

Step 4: Unplug the battery from its connector. Remove the battery and add the replacement battery at the location shown (see above picture) and position the wiring similar to the routing shown above. Plug the battery into its mating connector and reverse the steps noted above to complete the reassembly.

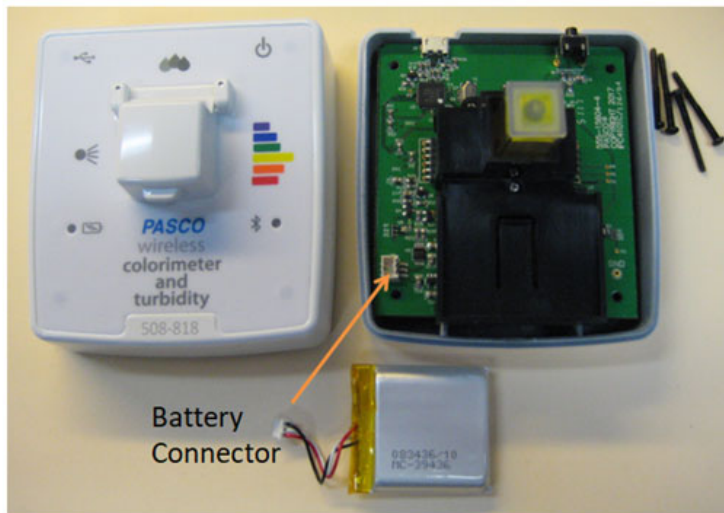

Battery Replacement and Disposal Instructions (Reference):

Batteries contain chemicals that, if released, may affect the environment and human health. Batteries should be collected separately for recycling, and recycled at a local hazardous material disposal location adhering to your country and local government regulations. To find a battery recycling drop-off location, please contact local waste disposal service or the product representative.



The battery or batteries used in PASCO products are marked with the European Union symbol for waste batteries (above) to indicate the need for separate collection and recycling. For small batteries, the symbol is printed on the packaging.

If you need further assistance, please contact Technical Support at support@pasco.com or through www.pasco.com on chat. To access chat through your browser, please enable all cookies on our website.

PASCO Form Number	Description	Date		Revision
013-16110	PS-3215 Battery Replacement Instructions	05-05-22	Page 3 of 3	B