

PS-3227; Wireless Sound Sensor Battery Replacement Instructions

Reference: Battery part number 540-113; Li-Poly 1000 mAH

Recommended Equipment: (PS-3299; Battery Replacement Tool Set)




Torx T5 Driver



Replacement Battery



CAUTION: Failure to use proper ESD protection could result in damage to the electronics.

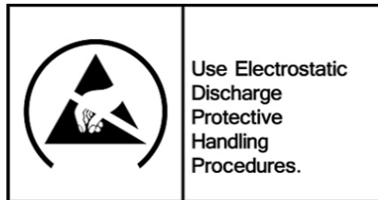

Step 1: Position the Sensor on a surface as shown. Note the location of the (4) Torx T5 Screws.



Torx T5 Screw, 4 Places

Step 2: Unscrew and remove the 4 Screws with the Torx T5 Screwdriver.



PASCO Form Number	Description	Date		Revision
013-17364	PS-3227 Battery Replacement Instructions	05-05-22	Page 1 of 3	A

Step 3

Separate the two halves and locate as shown.



Step 4:

Remove the battery hold down cover by removing the two screws with a Torx T5 Screwdriver. Unplug the battery and set aside.



PASCO Form Number	Description	Date	Page	Revision
013-17364	PS-3227 Battery Replacement Instructions	05-05-22	Page 2 of 3	A

Step 5 Place the new battery in the battery compartment and secure by reattaching the battery hold down cover. Plug in the battery into the battery connector and reverse the steps noted above to complete the reassembly.

Battery Replacement and Disposal Instructions (Reference):



Batteries contain chemicals that, if released, may affect the environment and human health. Batteries should be collected separately for recycling, and recycled at a local hazardous material disposal location adhering to your country and local government regulations. To find a battery recycling drop-off location, please contact local waste disposal service or the product representative.



The battery or batteries used in PASCO products are marked with the European Union symbol for waste batteries (above) to indicate the need for separate collection and recycling. For small batteries, the symbol is printed on the packaging.

If you need further assistance, please contact Technical Support at support@pasco.com or through www.pasco.com on chat. To access chat through your browser, please enable all cookies on our website.

PASCO Form Number	Description	Date		Revision
013-17364	PS-3227 Battery Replacement Instructions	05-05-22	Page 3 of 3	A