

PS-3232; Wireless //control.Node Battery Replacement Instructions

Battery part number: PS-3293; Li-Poly 3000 mAh

Recommended Equipment: (PS-3299; Battery Replacement Tool Set)




Torx T5 Driver



Replacement Battery



CAUTION: Failure to use proper ESD protection could result in damage to the electronics.

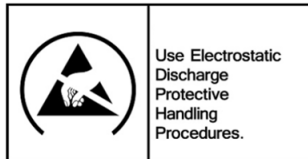

Step 1: Position the //control.Node looking at the bottom surface as shown. Note the location of the (4) Torx T5 Screws. Unscrew the (4) T5 Screws with a Torx T5 Screwdriver.



Torx T5 Screw, 4 Places

PASCO Form Number	Description	Date		Revision
013-17711	Battery Replacement Instructions	1-23-24	Page 1 of 4	A


Step 2: While holding the two halves together, flip the unit over looking at the top side as shown.



Step 3: Lift off the top cover and rotate it to the right as shown below. Gather the (4) screws and set aside. Note the orientation of the plug.



Step 4: Unplug the battery connector and locate the top cover as shown below. Unscrew the hold-down bracket T5 Screws (4 total) with a Torx T5 Screwdriver. Set the brackets and screws aside.

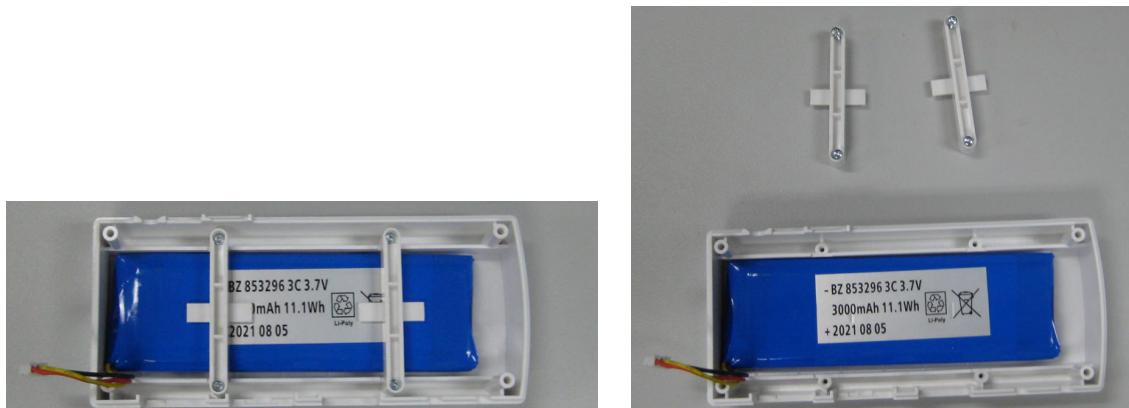


PASCO Form Number	Description	Date	Page 2 of 4	Revision
013-17711	Battery Replacement Instructions	1-23-24	Page 2 of 4	A

Step 5: Remove the old battery and replace it with the new replacement battery. Although the battery can be reinstalled in the cable up or cable down position, PASCO has determined that the cable up position is ideal for ease of re-assembly as shown below.



Step 6: Reassembly

- Replace the battery hold down brackets using the (4) T5 screws and the Torx T5 Screwdriver.
- Plug the battery connector into the main board (press firmly in place).
- Position the top cover as shown below and rotate it into place (this prevents the cable wires from being pinched).

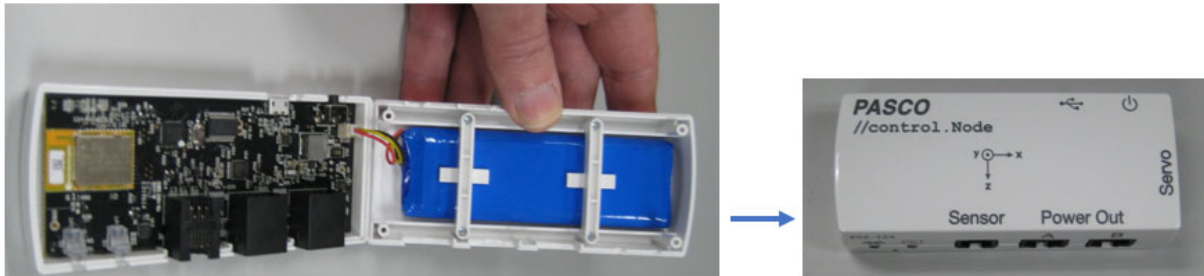

Step 7: Reassembly

While holding both halves together, flip the unit over and secure it with the (4) T5 screws and the Torx T5 Screwdriver. The Reassembly is complete.



Torx T5 Screw, 4 Places


PASCO Form Number	Description	Date		Revision
013-17711	Battery Replacement Instructions	1-23-24	Page 3 of 4	A

Battery Replacement and Disposal Instructions (Reference):

Batteries contain chemicals that, if released, may affect the environment and human health. Batteries should be collected separately for recycling and recycled at a local hazardous material disposal location adhering to your country and local government regulations. To find a battery recycling drop-off location, please contact local waste disposal service or the product representative.



The battery or batteries used in PASCO products are marked with the European Union symbol for waste batteries that indicate the need for separate collection and recycling. For small batteries, the symbol is printed on the packaging.

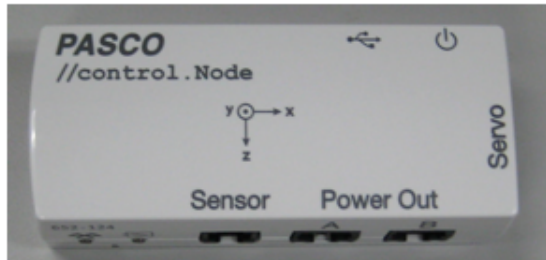

PASCO Form Number	Description	Date		Revision
013-17711	Battery Replacement Instructions	1-23-24	Page 4 of 4	A